

Phoenix 8000

User Manual

DOHRMANN™

Copyright

© 1997 by Tekmar-Dohrmann

No part of this booklet may be copied, reproduced, transmitted in any form or by any means, electronic or mechanical, including photocopying, recording, or information storage and retrieval systems, for any purpose other than the purchaser's personal use, without the express written permission of Tekmar-Dohrmann.

Updates

Tekmar-Dohrmann may update the information contained in this booklet without notice to purchaser.

The following are registered items:

Teflon®

Microsoft® Windows®

Microsoft® Excel®

DuPont® Corporation

Microsoft® Corporation

Microsoft® Corporation

Tekmar-DOHRMANN™

7143 East Kemper Road, Cincinnati, Ohio 45242-9576

Sales (800) 543-4461 • Outside the US and Canada (513) 247-7000 • Service (800) 874-2004

Telefax (513) 247-7050

Document Part Number: 14-7045-074

S:\Phoenix\Manual *.* .pm6 • V.02.12.97

Table of Contents

Preface

Understanding Safety Warnings	i
Working Safely	ii
Introduction to TOC Analysis	v
Analyzing TOC	vii

1 Introduction

Description of the Phoenix 8000	1-1
Unpacking the Phoenix 8000	1-3
Technical Specifications	1-4

2 Parts Description

Overview	2-1
Syringe Pump	2-2
IC Sparger	2-2
UV Reactor	2-3
Chlorine Scrubber	2-3
Moisture Control System	2-4
Nondispersive Infrared Detector (NDIR)	2-6
Flow Restrictors	2-7
Standard Flowmeters	2-7
Optional Flowmeters	2-7
Optional STS 8000 Autosampler	2-7

3 Preparing Reagents and Making Connections

Overview	3-1
Making Electrical Connections	3-1
Preparing Reagents	3-2
Understanding Gas, Water, and Reagent Connections	3-3
Tools and Supplies Needed for Connections	3-3
Swaging a Nut and Ferrule onto Tubing	3-4
Connecting Phoenix 8000 to a Gas Supply	3-5
Connecting Phoenix 8000 to a Water Supply	3-6
Connecting Phoenix 8000 to Acid and Persulfate Supplies	3-6
Filling Gas/Liquid Separator	3-7
Connecting Phoenix 8000 to Drain Line	3-7

Table of Contents

4 Connecting the Autosampler

Overview	4-1
Tools and Supplies Needed	4-2
Installing Fuses	4-2
Connecting RS-232 Cable	4-2
Setting Baud Rate	4-3
Removing Arm Locking Screw	4-3
Installing the Vertical Arm	4-4
Installing the Needle	4-5
Installing the Rinse Station and Support Bar	4-6
Installing the Tray	4-7
Installing the Rack	4-7
Connecting Power Supply	4-7

5 Understanding Operating Modes

Overview	5-1
Sample Introduction	5-1
Total Organic Carbon (TOC)	5-2
Total Carbon (TC)	5-5
Inorganic Carbon (IC)	5-7
Total Carbon Minus Inorganic Carbon (TC-IC)	5-9

6 Preparing to Analyze Samples

Overview	6-1
Installing the TOC Talk Software	6-1
Setting Flow Rates and Pressure	6-2
Understanding Calibration	6-3
Using TOC Talk	6-6
Instrument Setup	6-8
Diagnostics-Valves	6-9
Diagnostics-Syringe	6-10
Diagnostics-AutoSampler	6-12
Diagnostics-Communications	6-13
Diagnostics-Flowmeter Calibration	6-14
Calibration-Standards	6-15
Calibration-Set Active	6-17
Cal. Verification	6-18

(Chapter 6 - Preparing to Analyze Samples, cont.)

About 6-19
 Run Screen Choices 6-19
 Analysis Setup 6-27
 Method Setup Description of Terms 6-30
 Results 6-39
 Opening TOC Talk Files Using Microsoft Excel 6-48


7 Maintenance and Troubleshooting

Overview 7-1
 Daily Maintenance Checks 7-1
 Weekly Maintenance Checks 7-1
 Monthly Maintenance Checks 7-1
 Preventative Maintenance Check List 7-2
 Exterior Leak Checking of the Phoenix 8000 7-3
 Interior Leak Checking of the Phoenix 8000 7-5
 Calling Tekmar-Dohrmann Service 7-9
 Returning the Phoenix 8000 7-9
 Spare Parts List 7-10

Flow Diagram

Glossary

Index

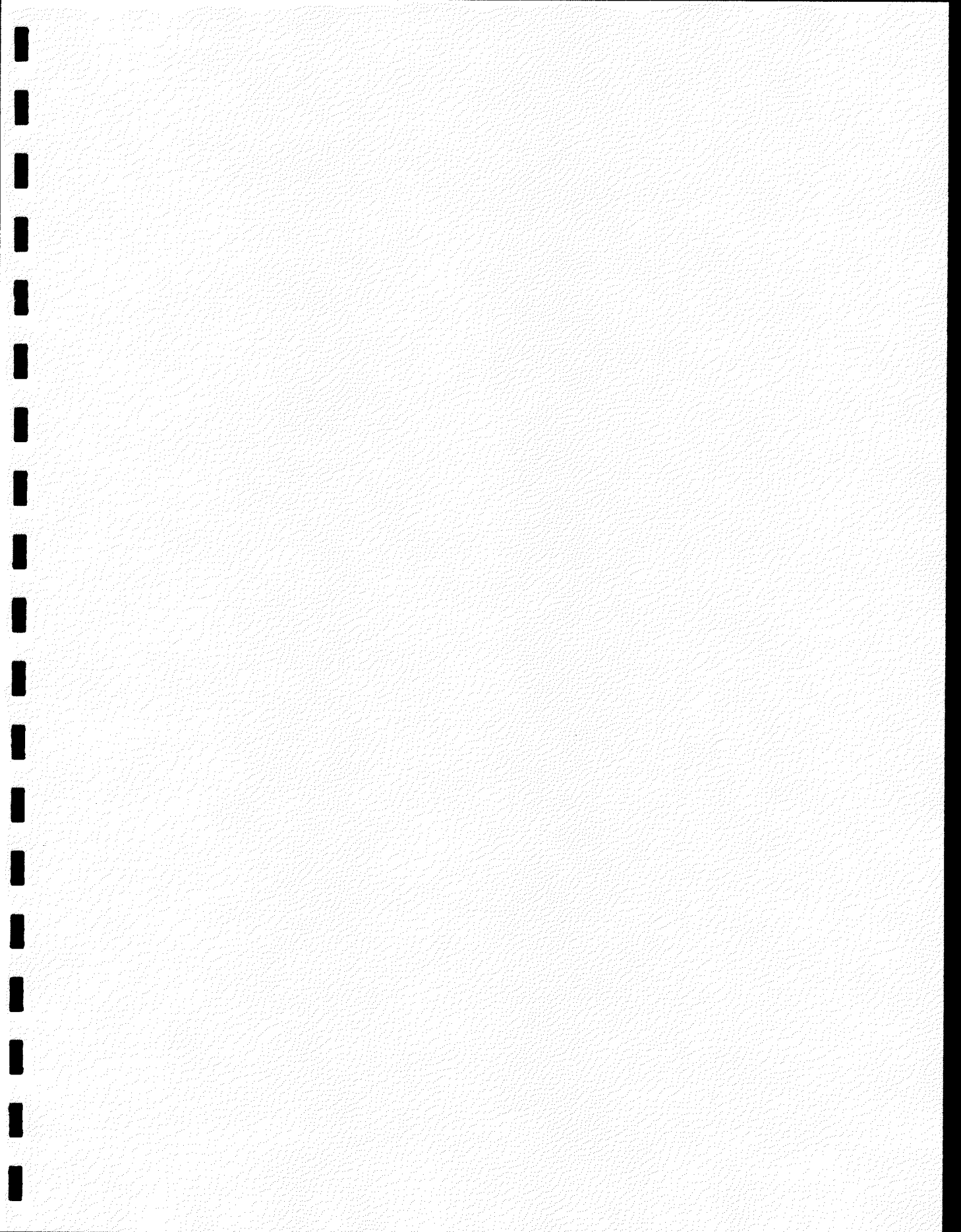

Understanding Safety Warnings

The table below describes symbols used in this manual.








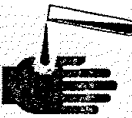




This symbol and/or notice	indicates:
 DANGER	imminently hazardous situation, which, if not avoided, will result in death or serious injury
 WARNING	potentially hazardous situation, which, if not avoided, can result in death or serious injury
 CAUTION	hazardous situation, which, if not avoided, will result in product or property damage and possible injury
	explosion
	hazardous voltage
	fire
	harmful fumes
	corrosive liquid or substance
	high temperature
	risk of eye injury; wear safety glasses
	hazardous or toxic substance
	moving parts can pinch and grab; keep hands and clothing away while power is on

Table 1 Explanation of Symbols