

Atomx Software Enhancement Installation

INSTRUCTIONS

1. If needed, backup all data prior to performing installation of software update. All old methods, custom methods, schedules, reports and calibrations will not be accessible after update is complete.
2. Uninstall the previous Atomx 1.0 software through the **Control Panel**, then **Add or Remove Programs**. (Figures 1A, 1B and 1C)
3. Download the Teklink Atomx file from the Extranet and place it in **C:\Program Files\Teledyne Tekmar**, or Insert the CD provided. If downloading from the Extranet, go to **C:\Program Files\Teledyne Tekmar\InstallMenu**. If using the installation CD, it should boot up automatically after a few moments. If not see your IT personnel.
4. Locate the **Atomx TekLink 1.1** file from the download or on the CD, left click the **Install** button and follow the on-screen prompts. (Figure 2)
5. Logon to the Atomx software.
6. Create new users and add new instruments as needed.



Figure 1A: Control Panel

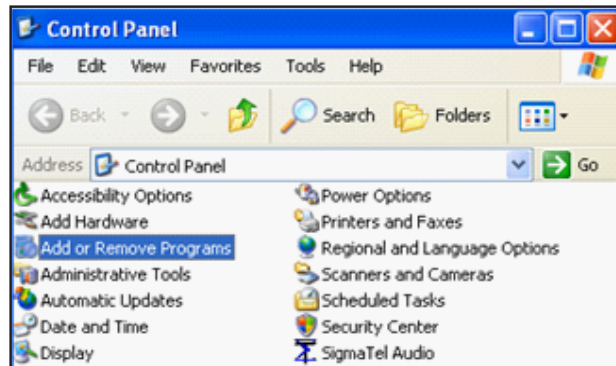


Figure 1B: Add or Remove Programs selected

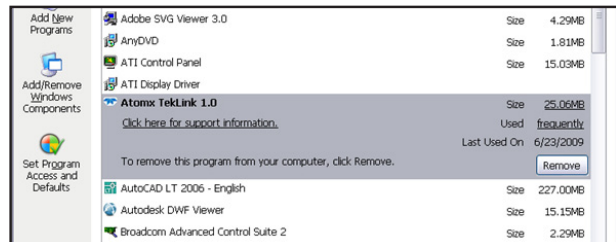


Figure 1C: Removal of Atomx Software

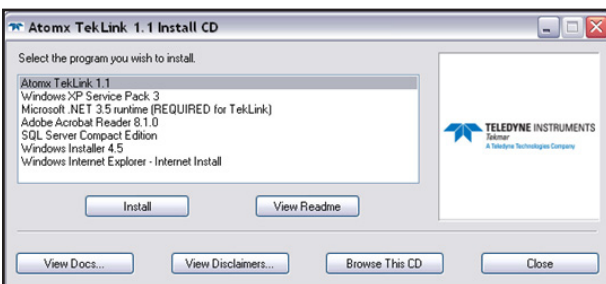


Figure 2: Installation of Atomx Teklink 1.1

Atomx Software Enhancement Installation

INSTRUCTIONS CONT'D

- Flash the firmware by going to **Tools** then **Update Firmware**. (Figure 3) Select **Browse**. If you downloaded the software, locate the file 102_1289.atomxmain.s19 from the Firmware folder on your C: drive. If using the CD, locate the file 102_1289.atomxmain.s19 from the Firmware folder and click **Open**.
- Once the appropriate file has been selected, click **Flash** and follow the on-screen prompts. Do not turn off the instrument during this time. (Figure 4)
- Once the Flashing of the firmware has been completed, power OFF the unit for 5 seconds then power ON to complete the installation.

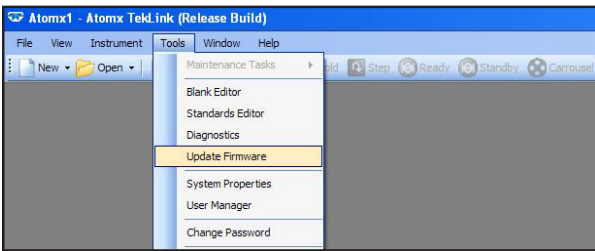


Figure 3: How to update through Tools | Firmware menu

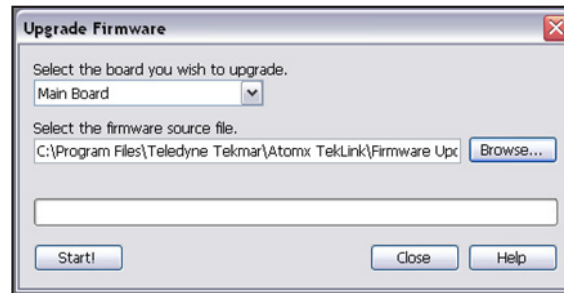


Figure 4: Click **Flash** to update the firmware and follow the on-screen prompts