

Installing the turret

- 1 For a 5890/6890 GC with a tray:
 - If you are changing from the 1-sample transfer turret to the 8-sample turret, disable the tray.
 - If you are changing from the 8-sample turret to the 1-sample transfer turret, enable the tray.

The 5890 Series II GC uses a software program run on a PC connected to the controller to configure the tray. Run the program, select **Configuration**, and either *check* (to enable) or *uncheck* (to disable) the tray. Select **Send Settings to ALS** to download the new setting to the controller.

All 6890 GCs disable the tray using the GC keypad.

- 2 Open the injector door. Unscrew and remove the knurled nut from the top of the turret.